



Syrris Ltd.
27 Jarman Way, Royston, Hertfordshire
SG8 5HW, United Kingdom
T: +44 (0)1763 242555
F: +44 (0)1763 242992
E: info@syrris.com
W: www.syrris.com

Syrris Limited
Registered office as address
Company No. 4257809

System/Product	Atlas
Model/Module	All
Subject Type	Baseware
Subject Version	All
Author	SV
Date	03/01/2008

Technical Note – How to check the version of Atlas base software (Baseware)

- a) To check the update has worked, select 'M' on the idle screen and press the atlas dial



- b) Select 'Administration'.



- c) Select 'Detected Hardware'.



- d) Select the 'Atlas Base'



- h) And finally, scroll down and check the atlas base version (Item 5 in the list)



Make a note of this version and let Support@Syrris.com know when requested.